



VINVENTIONS

WQS Color

Holder

Quick Start Guide

www.eu-vinventions.com

WQS Color

This document is intended to familiarize you with the NomaSense Color P100 holder.

Table of Contents

| | |
|-----------------------|---|
| 1. Packaging Contents | 2 |
| 2. Getting Started | 3 |
| 3. Maintenance | 4 |

1. Packaging Contents

| | |
|--------------------------|---------------------|
| 1. P100 Holder | 2. P100 Holder Cap |
| 3. 2x 5mm Cells | 4. 5mm Cell Adaptor |
| 5. Protective Carry Case | |



WQS Color

2. Getting Started




1. Remove the P100 Holder (1) and cap (2) from the case. Remove the cap (2).



2. Insert the 5 mm Cell Adaptor (4)



3. Turn on the P100, and calibrate as per the on screen instructions. Place into the P100 holder.

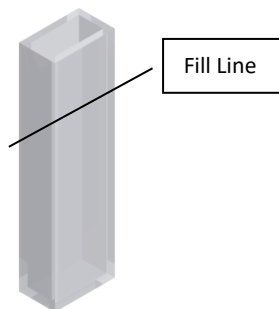

3. Fill the cell (3) with the required sample. Cell should be $\frac{2}{3}$ full.

WQS Color



4. Carefully slide the cell (3) into the P100 holder



5. Place on the P100 Holder Cap (2), and follow the P100 instructions to take a measurement.

3. Maintenance

Any spillage on to the white holder should immediately be wiped clean. Please use warm water with standard washing-up liquid and a soft cleaning cloth to achieve this. Do not use abrasive powders or harsh cleaning material on the holder.

Vinventions Sprl
ZI Les Plénesses
Chemin de Xhénorie, 7
4890 Thimister-Clermont
Belgium